


NOW AVAILABLE WITH  
**ANTI-FRINGING**  
BACK THINNED  
TECHNOLOGY

Spectroscopy

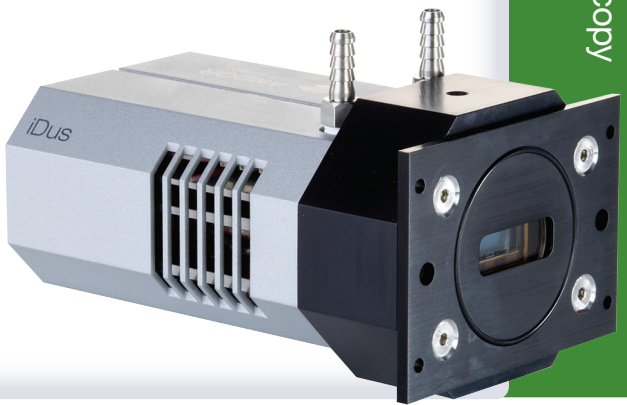

## Features and Benefits

- **Peak QE up to 95%**  
Visible-optimized 'BVF' & infrared-optimized 'BR-DD' model
- **Fringe suppression technology as standard**  
Fringing greatly eliminated (Deep-Depletion) or minimized (Back-Thinned 'BVF')
- **TE cooling down to -100°C**  
Critical for elimination of dark current detection limit - no inconvenience associated with LN<sub>2</sub>
- **Ultravac™ \*1**  
Critical for sustained vacuum integrity and to maintain unequalled cooling and QE performance, year after year
- **Single UV-grade fused silica window**  
Best UV-NIR throughput performance, specific AR coating and wedge options available
- **26 x 26 µm pixels**  
Optimized format for high dynamic range and resolution
- **Software-selectable pre-amplifier gain**  
Choice of best SNR performance or dynamic range at the touch of a button
- **USB 2.0 connection**  
Ideal for laptop operation  
Seamless operation alongside USB-based Shamrock spectrograph family
- **Compact and rugged platform**  
Ideal for OEM integration & laboratory setups
- **Solis software for Spectroscopy**  
Comprehensive, user-friendly interface for simultaneous detector & spectrograph control
- **Software Development Kit (SDK)**  
Ease of control integration into complex setups: Matlab, Labview, Visual Basic or C/C++

## Workhorse laboratory and OEM CCD platform

The USB 2.0-based iDus series is a compact, yet feature-rich platform suitable for demanding spectroscopy applications such as low-light UV/NIR Photoluminescence or Raman spectroscopy, as well as day-to-day routine laboratory operation and integration into industry-grade systems.

The iDus 401 boasts market-leading deep thermo-electric cooling to -100°C, complemented by Andor's Ultravac™ technology with its unmatched reliability track record in the scientific and industrial community. The iDus low-noise electronics allows seamless access to the best signal-to-noise ratio performance for the widest range of photon detection regimes.

**Anti-fringing comes as standard** on back-illuminated sensors for this 1024 x 127, 26 µm pixels array optimized for high dynamic range and high resolution acquisitions from far UV to NIR.

## Specifications Summary \*2

Active pixels	1024 x 127
Pixel size (W x H)	26 x 26 µm
Image area	26.6 x 3.3 mm
Register well depth (typical)	1,000,000 e <sup>-</sup>
Maximum cooling	-100°C
Maximum spectra per sec	88
Read noise	As low as 3 e <sup>-</sup>
Dark current	As low as 0.0005 e <sup>-</sup> /pixel/sec

## Key Specifications \*2

Model number	DV401A	DU401A	DU401A BR-DD
Sensor options	<ul style="list-style-type: none"> <li>• BVF: Back Illuminated CCD, Vis-optimized and anti-fringing</li> <li>• FI: Front Illuminated CCD</li> </ul>	<ul style="list-style-type: none"> <li>• BVF: Back Illuminated CCD, Vis-optimized and anti-fringing</li> <li>• FI: Front Illuminated CCD</li> </ul>	<ul style="list-style-type: none"> <li>• BR-DD: Back Illuminated CCD, Deep Depletion with anti-fringing</li> <li>Optimum sensor for Near IR applications</li> </ul>
Active pixels *3	1024 x 127		
Pixel size	26 x 26 $\mu\text{m}$		
Image area	26.6 x 3.3 mm with 100% fill factor		
Minimum temperatures *4			
Air cooled	-55°C	-80°C	-80°C
Coolant recirculator	-65°C	-95°C	-95°C
Coolant chiller, coolant @ 10°C, 0.75l/min	-70°C	-100°C	-100°C
Max spectra per second *5	88 (10 row crop mode), 81 (Full Vertical Bin)		
System window type	Single UV-grade fused silica window, uncoated. Various AR coatings & MgF <sub>2</sub> options available	Single UV-grade fused silica window, uncoated. Various AR coatings & MgF <sub>2</sub> options available	Single UV-grade fused silica window; 1° wedge, AR coated on both sides, optimized at 900 nm
Blemish specifications	Grade 1 as per sensor manufacturer definition		

## Advanced Specifications \*2

Dark current, e <sup>-</sup> /pixel/sec @ max cooling									
FI	0.0035			0.0005			-		
BVF	0.006			0.003			-		
BR-DD	-			-			0.013		
Register well depth	1,000,000 e <sup>-</sup>								
Read noise (e <sup>-</sup> ) *6	33 kHz	50 kHz	100 kHz	33 kHz	50 kHz	100 kHz	33 kHz	50 kHz	100 kHz
FI: Typ (Max)	3 (5)	4 (6)	8 (12)	3 (5)	4 (6)	9 (12)	-	-	-
BVF: Typ (Max)	7 (8)	10 (12)	12 (15)	6 (8)	10 (12)	12 (15)	-	-	-
BR-DD: Typ (Max)	-	-	-	-	-	-	5 (8)	6 (12)	12 (15)
Sensitivity (e <sup>-</sup> /count)	33 kHz	50 kHz	100 kHz	33 kHz	50 kHz	100 kHz	33 kHz	50 kHz	100 kHz
High Dynamic Range mode	2	3.5	14	2.5	3.5	16	2.5	3.5	20
High Sensitivity mode	-	2.5	9	-	2.5	11	-	2.5	12
Linearity *7	Better than 99%								
Digitization	16 bit								
Vertical clock speed *8	8, 16, 32, 64 $\mu\text{s}$ (software selectable)								

## Have you found what you are looking for?

**Need a larger sensor?** The iDus 420 series is similar to the 401 series with 26.6 x 6.6 mm format and 1024 x 255 pixels.

**Need to work further into the NIR?** The iDus InGaAs series, with up to 1024 pixel linear array with transmission to 2.2  $\mu\text{m}$ .


**Need a faster or more sensitive CCD?** The Newton series offers over 1,600 spectra per second and EM technology.

**Need more UV sensitivity?** The iDus 420 series is similar to the 401 series with a range of UV optimized sensors.


**Need a customized version?** Please contact us to discuss our Customer Special Request options.

The iDus series combines seamlessly with Andor's research grade Shamrock Czerny-Turner spectrographs.


## Dark Current <sup>\*9</sup>

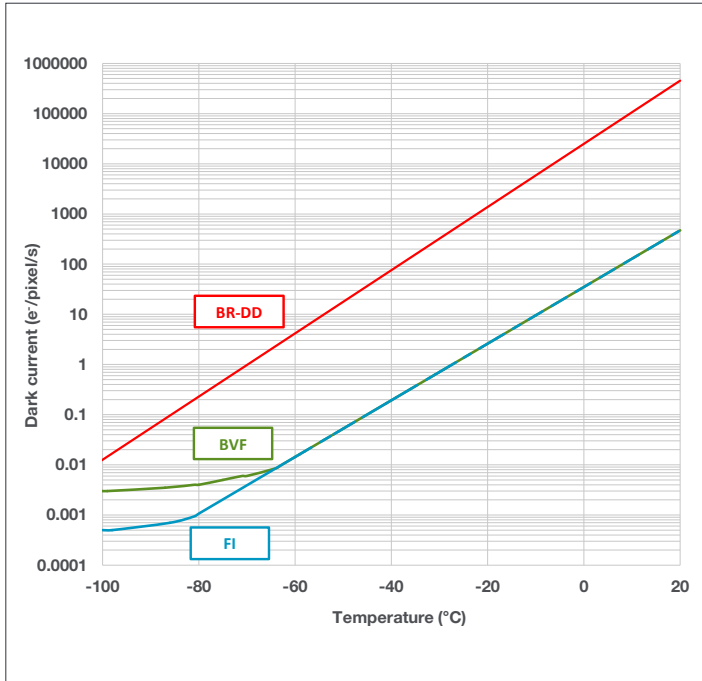


## Quantum Efficiency Curves <sup>\*10</sup> 25°C

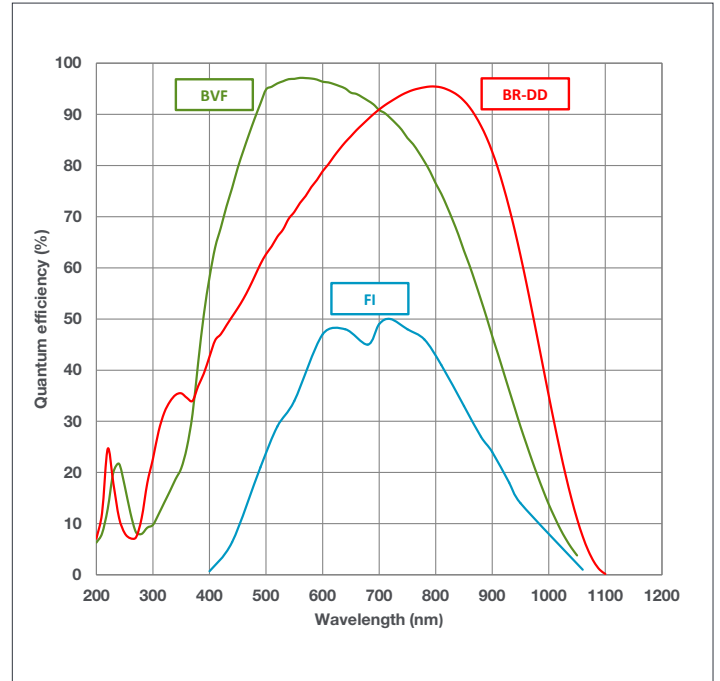

## Typical Setup

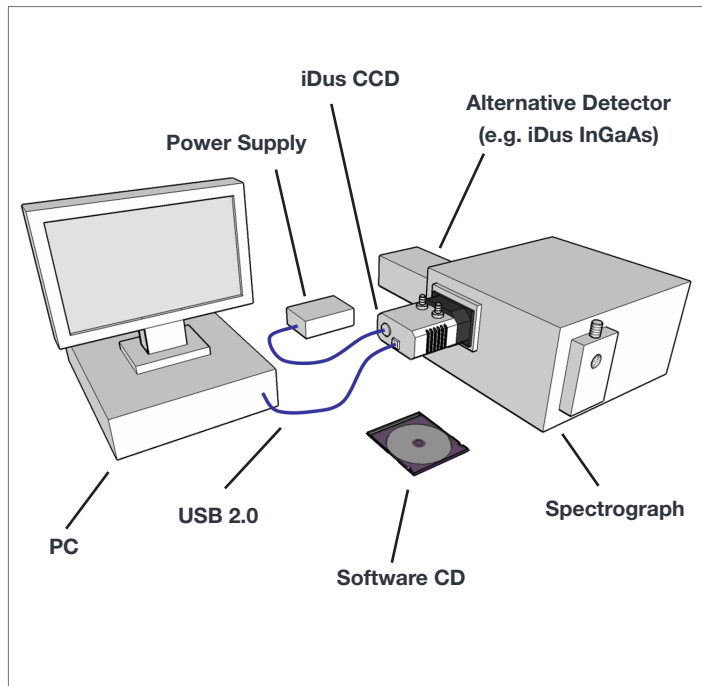

## Quantum Efficiency Curves <sup>\*10</sup> -100°C

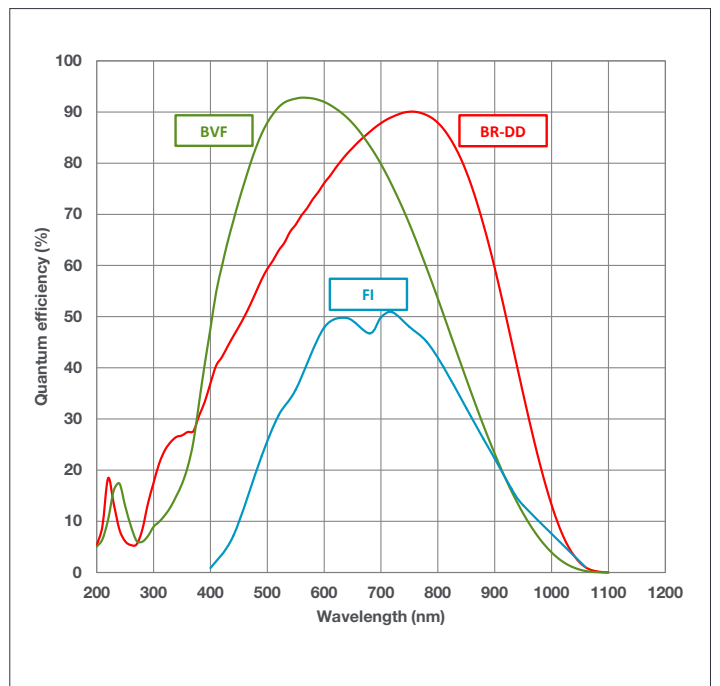

## Creating The Optimum Product for You

How to customize the iDus 401A:

### Step 1.

Simply select from the 2 cooler types that best suit your needs from the selection opposite.

### Step 2.

The iDus 401A comes with 3 options for sensor types. Please select the sensor which best suits your needs.

### Step 3.

Please indicate if you wish to select an alternative window and which software you require.

### Step 4.

For compatibility, please indicate which accessories are required.



iDus CCD mounted on a Shamrock 500 mm triple grating imaging spectrograph, ideal for high resolution spectroscopy.



### Step 1.

#### Choose cooler type

**DV:** Standard cooling, to -70°C (not a valid option with the BR-DD model)  
**DU:** Deep cooling, to -100°C

### Step 2.

#### Choose sensor type

**BR-DD:** Back Illuminated, Deep Depletion CCD with anti-fringing  
**BVF:** Back Illuminated CCD, Vis-optimized with anti-fringing  
**FI:** Front Illuminated CCD

### Step 3.

**The 401 FI & BV models are supplied with an uncoated UV-grade fused silica window as standard. BR-DD models are provided with an NIR optimized AR coating. The following alternative window choices are available and must be ordered at time of build (if selected):**  
**OPTION-C1-AR1** AR coated UV-grade fused silica window (optimized broadband visible 400-900nm). 50% transmission at 180 nm  
**OPTION-C1-MGF2** Magnesium Fluoride window for transmission in the VUV. 50% transmission at 120 nm

#### The iDus CCD also requires at least one of the following software options:

**Solis for Spectroscopy** A 32-bit application compatible with 32 and 64-bit Windows (XP, Vista, 7 and 8) offering rich functionality for data acquisition and processing. AndorBasic provides macro language control of data acquisition, processing, display and export. Control of Andor Shamrock spectrographs and a very wide range of 3<sup>rd</sup> party spectrographs is also available, see list below.

**Andor SDK** A software development kit that allows you to control the Andor range of cameras from your own application. Available as 32 and 64-bit libraries for Windows (XP, Vista, 7 and 8) and Linux. Compatible with C/C++, C#, Delphi, VB6, VB.NET, LabVIEW and Matlab.

### Step 4.

#### The following accessories are available:




**XW-RECR** Coolant re-circulator for enhanced cooling performance  
**ACC-XW-CHIL-160** Oasis 160 Ultra Compact Chiller Unit (tubing to be ordered separately)  
**ACC-6MM-TUBING-2xxxxM** 6 mm tubing option for ACC-XW-CHIL-160  
**LM-C** C-mount lens adaptor  
**LM-NIKON-F** C-mount lens adaptor  
**LMS-NIKON-F-NS25B** Nikon F-mount lens adaptor with shutter  
**ACC-SD-VDM1000** Shutter Driver for NS25B Bistable Shutter (not needed for Shamrock spectrographs)  
**ACC-SHT-NS25B** Bistable Shutter, Standalone (not needed for Shamrock spectrographs)

#### Spectrograph Compatibility


The iDus series is fully compatible with Andor's Shamrock spectrograph (163 - 750 nm focal lengths) family. Spectrograph mounting flanges and software control are available for a wide variety of 3<sup>rd</sup> party spectrographs including McPherson (including 1 m and greater focal length option), JY/Horiba (excluding USB models), PI/Acton, Chromex/Bruker, Oriel/Newport, Photon Design, Dongwoo, Bentham, Solar TII and others.

## Product Drawings


Dimensions in mm [inches]

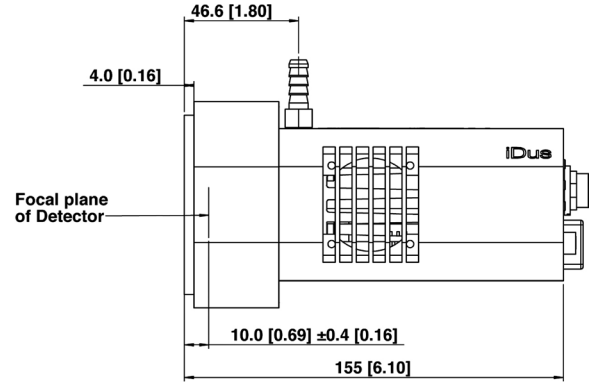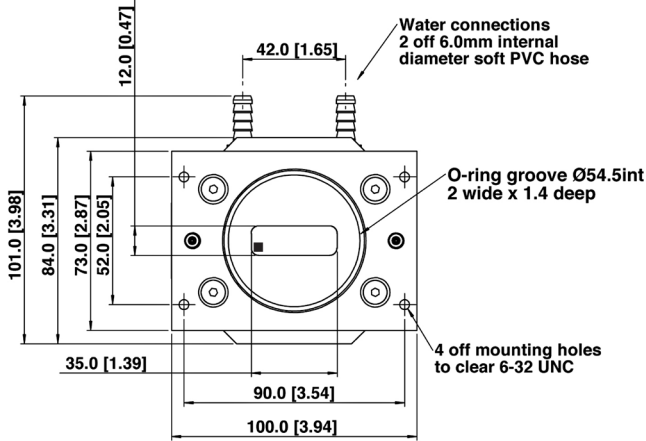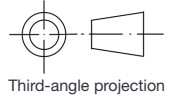

■ = position of pixel 1,1  
Weight: 2 kg [4 lb 8 oz]

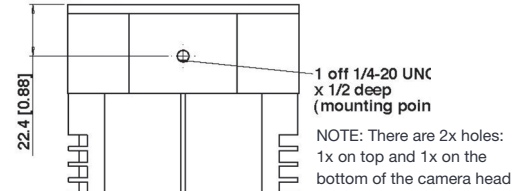

Mounting hole locations

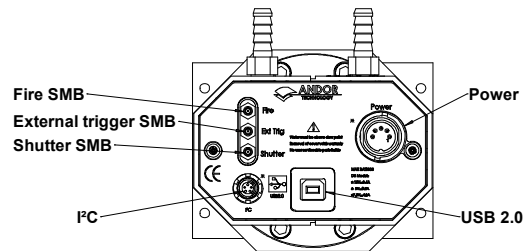

Rear connector panel

## Connecting to the iDus

### Camera Control

Connector type: USB 2.0

### TTL / Logic

Connector type: SMB, provided with SMB - BNC cable  
1 = Fire (Output), 2 = External Trigger (Input), 3 = Shutter (Output)

### I<sup>2</sup>C connector

Compatible with Fischer SC102A054-130  
1 = Shutter (TTL), 2 = I<sup>2</sup>C Clock, 3 = I<sup>2</sup>C Data, 4 = +5 Vdc, 5 = Ground

### Minimum cable clearance required at rear of camera

90 mm

## Applications Guide

	DV 401 FI	DV 401 BVF	DU 401 FI	DU 401 BVF	DU 401 BR-DD
Absorption/Transmittance/Reflection	✓	✓	✓	✓	✓
Atomic Emission Spectroscopy				✓	
Fluorescence & Luminescence	✓	✓	✓	✓	✓
NIR Spectroscopy			✓		✓
Raman Spectroscopy (244 – 488 nm)	✓	✓		✓	
Raman Spectroscopy (514, 532, 633 nm)	✓	✓	✓	✓	
Raman Spectroscopy (785, 830 nm)	✓		✓		✓

✓ = Suitable  
✓ = Optimum



## Order Today

Need more information? At Andor we are committed to finding the correct solution for you. With a dedicated team of technical advisors, we are able to offer you one-to-one guidance and technical support on all Andor products. For a full listing of our local sales offices, please see: [andor.com/contact](http://andor.com/contact)

### Our regional headquarters are:

#### Europe

Belfast, Northern Ireland  
Phone +44 (28) 9023 7126  
Fax +44 (28) 9031 0792

#### Japan

Tokyo  
Phone +81 (3) 3518 6488  
Fax +81 (3) 3518 6489

#### North America

Connecticut, USA  
Phone +1 (860) 290 9211  
Fax +1 (860) 290 9566

#### China

Beijing  
Phone +86 (10) 5129 4977  
Fax +86 (10) 6445 5401

### Items shipped with your camera:

- 1x 2m BNC - SMB connection cable
- 1x 3m USB 2.0 cable Type A to Type B
- 1x Set of Allen keys (7/64" & 3/32")
- 1x Power supply with mains cable
- 1x Quick launch guide
- 1x CD containing Andor user guides
- 1x Individual system performance booklet
- 1x CD containing either Solis software or SDK (if ordered)

### Footnotes: Specifications are subject to change without notice

1. Assembled in a state-of-the-art facility, Andor's UltraVac™ vacuum process combines a permanent hermetic vacuum seal (no o-rings), with a stringent protocol and proprietary materials to minimize outgassing. Outgassing is the release of trapped gases that would otherwise degrade cooling performance and potentially cause sensor failure.
2. Figures are typical unless otherwise stated.
3. Edge pixels may exhibit a partial response.
4. Minimum temperatures listed are typical values with ambient temperature of 20°C.
5. Based on a Horizontal Pixel Readout of 100 kHz and a vertical pixel shift of 8 μs.
6. System Readout Noise is for the entire system. It is a combination of CCD readout noise and A/D noise. Measurement is for Single Pixel readout with the CCD at a temperature of -50°C (BR-DD models tested at -80°C) and minimum exposure time under dark conditions. Noise is measured at the highest available pre-amplifier gain for each speed.
7. Linearity is measured from a plot of counts vs exposure time under constant photon flux up to the saturation point of the system.
8. Vertical speeds are software selectable. All sensors are designed to give optimum ChargeTransfer Efficiency (CTE) at 16 μs vertical pixel shift, some decrease in CTE may be observed at faster shift speeds.
9. The graph shows typical dark current level as a function of temperature. The dark current measurement is averaged over the CCD area excluding any regions of blemishes.
10. Quantum efficiency of the sensor as supplied by the sensor manufacturer.

### Minimum Computer Requirements:

- 3.0 GHz single core or 2.4 GHz multi core processor
- 2 GB RAM
- 100 MB free hard disc to install software (at least 1 GB recommended for data spooling)
- USB 2.0 High Speed Host Controller capable of sustained rate of 40 MB/s
- Windows (XP, Vista, 7 and 8) or Linux

### Operating & Storage Conditions

- Operating Temperature: 0°C to 30°C ambient
- Relative Humidity: < 70% (non-condensing)
- Storage Temperature: -25°C to 50°C

### Power Requirements

- 110 - 240 VAC, 50 - 60 Hz



Windows is a registered trademark of Microsoft Corporation.  
Labview is a registered trademark of National Instruments.  
Matlab is a registered trademark of The MathWorks Inc.

SiDus401SS 0713 R1